LN₂ for OPEN TOP VESSELS SOP (NOT FOR DEWARS)



September 2020

NOTE: ALWAYS wear PPE when working with LN₂.

Before filling your vessel, you MUST do the following:

- **1.** Wear long pants.
- **2.** Have a buddy. No one is permitted to fill a vessel alone.
- 3. Scan your entry card for the number of vessels you are filling.
- **4.** Remain with the vessel to be filled at all times during the fill process.
- **5.** Wear the provided PPE before starting the fill process. PPE consists of a face shield, hearing protection and blue cryogenic gloves.

FILLING PROCEDURE

- 1. Make sure that the diffuser is on the end of flex line. If it is not there, contact MNTC staff immediately. **DO NOT FILL A VESSEL WITHOUT THE DIFFUSER.**
- 2. Place the flex line on the concrete away from your body.
- 3. Open the small valve on the fill line.
- **4.** Open the larger valve by the LN2 tank.
- 5. Wait for a flow of LN2 to exit from the diffuser to prime the line with LN2.
- 6. Turn off the small valve on the fill line.
- **7.** Place the flex into the top of your vessel.
- 8. Open the small valve on the fill line and fill the dewar to your desired level.
- **9.** When finished close the small valve on the fill line.
- **10.** Close the large valve on the LN2 tank, but **DO NOT OVERTIGHTEN**.
- 11. Place flex end on concrete and away from your body.
- **12.** Open the small valve on the fill line to purge remaining LN₂ in the line. Keep the line open until you hear a high pitched whine.
- 13. Close the small valve, but do not overtighten.
- **14.** Place the PPE back into the provided box and close the lid on the box.